



**NOTES:**

1. THE ANSI Z358.1-2009 STANDARD SPECIFIES EYEWASH NOZZLE HEIGHTS TO BE 33 IN. TO 45 IN. ABOVE THE FLOOR AND 6 IN. MINIMUM FROM WALL OR NEAREST OBSTRUCTION.
2. INSTALL SO THAT EYEWASH HEAD IS OVER SINK WHEN ACTIVATED. EYEWASH REQUIRES 4-1/2" OF CLEARANCE AROUND HEAD FOR FACEWASH STREAMS.
3. DRILL ONE Ø7/8" HOLE AND ONE Ø3/32" HOLE WHERE UNIT IS TO BE INSTALLED. REFER TO MOUNTING HOLE DETAIL. GO TO P102C OR THE LATEST REVISION FOR SHOWER DETAIL REFERENCE.
4. THESE DIMENSIONS MAY VARY ±1/2".
5. ALL SUBSTITUTE MUST BE APPROVED BY LBNL PROJECT MANAGER & EH&S.
6. SEE P102C OR THE LATEST REVISION FOR SHOWER DETAIL AND ELECTRICAL REQUIREMENT.

**SINGLE ACTION PULL DOWN MODEL 7610  
EMERGENCY EYE/FACEWASH & SHOWER INSTALLATION DETAIL**

SCALE: NONE

X  
XX

LBNL STANDARD  
EMERGENCY EYE/FACEWASH & SHOWER INSTALLATION DETAIL  
SINGLE ACTION PULL DOWN MODEL 7610

UNIVERSITY OF CALIFORNIA LAWRENCE BERKELEY NATIONAL LABORATORY  
FACILITIES DIVISION

DRAWN BY	RZT	DATE	12/22/2011
CHECKED BY	MCD		12/22/2011
APPROVED BY	MCD		12/22/2011
CAD FILE PATH			
SCALE	NTS		
DRAWING NO.	P110A		SHEET
PROJECT NO.			